



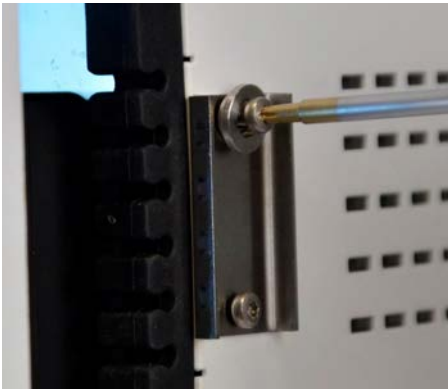
## Mounting bracket installation for AZURA® VU 4.1

### Product information

The intended use of the mounting bracket is to install the valve drive AZURA Valve Unifier VU 4.1 on the side of any AZURA L device. The following instruction depicts the installation at the right side of an AZURA L device but is applicable for the left side as well.

### Procedure

**Tools** Torx key, size TX10

Steps	Figure
1. Remove the screws on the left side of the valve drive.	
2. Align the mounting bracket with the bore holes of the valve drive and screw it tight with the provided torx screws.	
3. Align the guidance clamp with the bore holes of the right-side panel of the AZURA L device and screw it tight with the provided torx screws. Make sure that the upper screw is installed together with the washer.	

## Steps

- Put the upper keyhole of the mounting bracket onto the washer at the AZURA L device (1) and slide down the valve drive until the middle hole locks with the lower screw of the guidance clamp (2).

