Produktbeschreibung AZURA Click

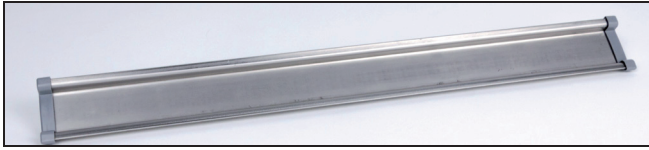
Abb. 1 Tragschiene AZURA Click

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AZURA Click

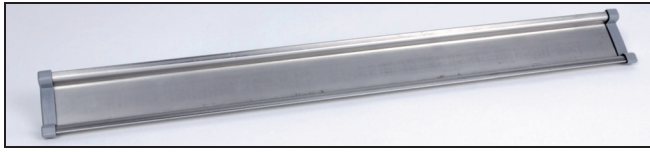
Fig. 1 Carrier rail AZURA Click

The carrier rail AZURA Click is for mounting an interface box, the Pressure Control or the Air Sensor. The carrier rail will be stuck to the side panel of your AZURA L device in 9 cm height.

Attaching

Precondition

- The tubular film is removed.
- The end caps are attached.
- The side panel is clean, dry, and free from dust and grease.

Tools

Measuring instrument with millimeterscale for linear measurement (e. g. folding rule, ruler o. ä.).

Note: Do not stick the carrier rail over the ventilation grille.

Process

1. Remove the protective foil (Fig.2).

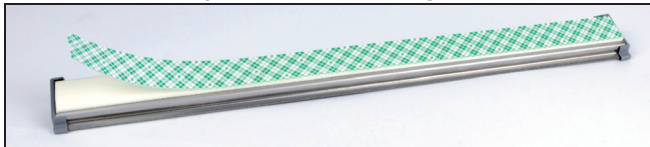
Fig. 2 Remove the protective foil

2. Position the carrier rail flush with the back panel in 9 cm height (Fig.3 alike).
3. Push the carrier rail strong against the side panel.

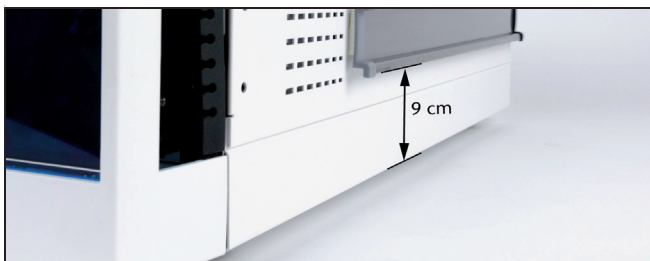
Fig. 3 Position of the carrier rail

Mounting the Holding Device at the Interface Box

Tools

Crosstip screwdriver size PH1.

Process

1. Position the holding device with the handle down at the interface box.
2. Screw the socket in the holding device and the interface box.

Attaching the Interface Box

Precondition

- The holding device is fixed.
- AZURA Click is attached.

Process

1. Hook the interface box on the upper border of the carrier rail (Fig.4).
2. Push the interface box against the lower border of the carrier rail.

Fig. 4 Position of the interface box

Taking off the Interface Box

Process

1. Pull the handle of the holding device down ① (Fig.5).
2. Take off the interface box sloping upwards ②.

Fig. 5 Take off the interface box

Attaching the Pressure Control

Precondition

- AZURA Click is attached.

Process

1. Hook the Pressure Control on the upper border of the carrier rail.
2. Push the Pressure Control against the lower border of the carrier rail.

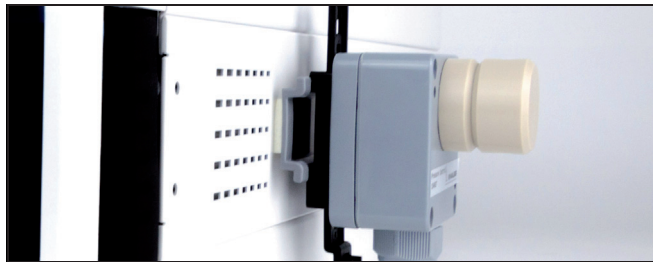
Fig. 6 Position of the Pressure Control

Taking off the Pressure Control

To take off the Pressure Control go ahead, like written before in "Take off the interface box".

Attaching of the Air Sensor

Precondition

- The metal rail is removed (Fig.7).
- AZURA Click is attached.

Process

1. Pull back the metal rail against the spring.
2. Take off the metal rail.

Fig. 7 Take off the metal rail

3. Push the Air Sensor sloping against the lower border of the carrier rail (Fig.8).
4. Pull the Air Sensor to the top.
5. Push the Air Sensor over the upper border of the carrier rail

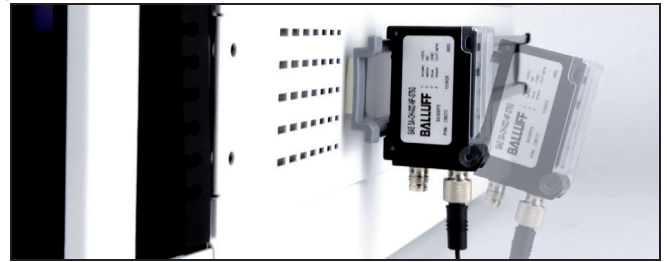
Fig. 8 Position of the Air Sensors

Taking off the Air Sensor

Process

1. Pull the Air Sensor over the top of the upper carrier rail.
2. Take off the Air Sensor.

Repeat Orders

Name	Order no.
AZURA Click	A70089
Interface Box	AZB00
Air Sensor for 1/16" Tubing	A70092-1
Air Sensor for 1/14" Tubing	A70093-1
Pressure Control	AZG10

Technical Data

Dimensions (B x H x T)	330 x 35 x 7,5 mm
Material	Stainless steel