




**Note:** Read the corresponding technical documentation for handling and safety reasons.

## Installation capillaries 1/8"

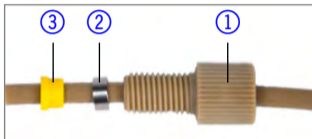

1. Ensure that the capillary ends are flat and even.
2. Push the capillary through the head of the nut ① .
3. Slide the seal ring ② over the capillary. The tapered end has to point away from the nut.
4. Attach the ferrule ③ .
5. Fasten the capillary hand-tight.



**KNAUER**  
**Wissenschaftliche Geräte GmbH**  
**Hegauer Weg 38**  
**14163 Berlin, Germany**

**Phone: +49 30 809727-0**

**Fax: +49 30 8015010**

**E-Mail: [info@knauer.net](mailto:info@knauer.net)**

**Internet: [www.knauer.net](http://www.knauer.net)**