




Read the corresponding technical documentation for handling and safety reasons.

## Installation




Ensure that the capillary ends are flat and even.

1. Push the capillary through the head of the nut ①.
2. Attach the ferrule ②.

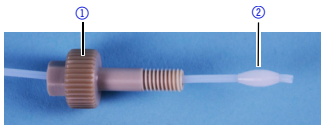

3. Fasten the capillary at the flow cell ③ fingertight.

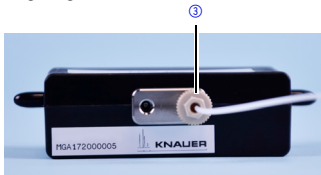



**KNAUER**

**Wissenschaftliche Geräte GmbH**

**Hegauer Weg 38**

**14163 Berlin, Germany**

**Phone: +49 30 809727-0**

**Fax: +49 30 8015010**

**E-Mail: [info@knauer.net](mailto:info@knauer.net)**

**Internet: [www.knauer.net](http://www.knauer.net)**